

Pre-made anti-microRNA (for human and mouse) lentivirus User Manual

Catalog#	Product Name		Amounts
A-miRNA H1-Bsd	anti miRNA (H1) with Blasticidin marker		
A-miRNA H1-GB	anti miRNA (H1) with GFP-Blasticidin marker	1	0.5 ml of outin
A-miRNA H1-GP	anti miRNA (H1) with GFP-Puromycin marker	1.	0.5 mi or anti a
A-miRNA H1-Puro	anti miRNA (H1) with Puromycin marker		microPNA
A-miRNA H1-RB	anti miRNA (H1) with RFP-Blasticidin marker		lentivirus
A-miRNA H1-RP	anti miRNA (H1) with RFP-Puromycin marker		icituviius,
A-miRNA U6-Bsd	Anti miRNA (U6) with Blasticidin marker	2	0.5ml of Negative
A-miRNA U6-GB	Anti miRNA (U6) with GFP-Blasticidin marker	2.	control anti-
A-miRNA U6-GP	Anti miRNA (U6) with GFP-Puromycin marker		microRNA
A-miRNA U6-Puro	Anti miRNA (U6) with Puromycin marker		lentivirus
A-miRNA U6-RB	Anti miRNA (U6) with RFP-Blasticidin marker		
A-miRNA U6-RP	Anti miRNA (U6) with RFP-Puromycin marker		

Storage: < -70 °C, avoid repeat freeze/thaw cycles. Stable for > 6 months.

Product Description:

1. Introduction:

MicroRNAs (**miRNAs**) are small non-coding RNA molecule found in plants and animals. They are transcribed from RNA precursors (**pre-miRNAs**), and have a mature length of 19-26 nucleotides. MiRNAs silence or repress gene expression by binding to complementary sequences within gene-coding mRNAs. Each miRNA can target multiple genes, and it is believed that >60% of mammalian gene expression is regulated by miRNAs. This includes most biological processes, including a variety of diseases and developmental disorders. More than 2000 mature miRNAs have been discovered in human and more than 1200 in mouse so far.

Anti-miRNAs, or miRNA inhibitors are short RNA molecules that are complementary to mature miRNAs. Anti-miRNA specifically binds to endogenous miRNA, thereby inhibiting or down-regulating miRNA activity. For this reason, anti-miRNAs are very useful tools for loss-of-function research involving specific miRNAs. Anti-miRNAs are also potential therapeutic agents for some diseases.





Anti-miRNA can be delivered by two methods: 1) single stranded nucleic acids designed to specifically bind the target miRNA sequence, or 2) vector-based expression of the antisense sequence.

GenTarget's ready-to-use, anti-miRNA lentiviruses are produced from our optimally designed anti-miRNA lentivectors. The annealed DNA oligonucleotide hairpin duplex encoding the antisense sequence of a specific miRNA was cloned under a **constitutive** human U6 promoter or under an <u>optional inducible</u> human H1 promoter. GenTarget's anti-miRNA lentivectors have been carefully designed for the highest anti-miRNA expression level, precise promoter transcription position, and favorability of the antisense strand for RISC complex processing. The engineered asymmetric duplex transcribes a perfectly matched antisense RNA molecule and a leftover bulged-out sense RNA molecule. The antisense RNA molecule is expressed at a high level and complementarily "binds" to and inactivates endogenous, mature miRNA. Please see the map below for the anti-miRNA lentivector core structure.



2. **Pre-made anti-miRNA lentivirus**:

Anti-miRNA lentivirus is the easiest and most reliable method for consistent delivery of anti-miRNA molecules. GenTarget's anti-miRNA lentiviruses feature **a choice of two promoters** (H1 or U6) and **a variety of selection markers** (antibiotic markers or fluorescent-antibiotic fusion markers), providing you with tools for long-term constitutive or inducible inhibition of targeted miRNA functionality. Lentivirus can be transduced into the majority of mammalian cell types, and the human H1 and U6 promoters are active in almost all mammalian cells. GenTarget provides packaged anti-miRNA lentivirus for any specific miRNA from human, mouse, or any other species.



Click here to see GenTarget's anti-human miRNA list and anti-mouse miRNA list.

All lentiviruses demonstrate strong transduction efficiency, and each lot of virus is validated for titer of approximately $1 \times 10^{7-8}$ IFU/ml. Quality is guaranteed.

Pre-made lentivirus is provided as ready-to-use **0.5ml** aliquots in DMEM medium with 10% FBS and 60μ g/ml polybrene. Upon request, lentivirus can be provided in PBS solution *in vivo* use.

For general questions about our ready-to-use lentiviral particles, please see **FAQ for pre-made lentiviral particles** (.pdf) on our website.

Note: GenTarget also provides pre-miRNA expression lentivirus for any specific miRNA; please contact us for a service quote.

3. What is the negative control for anti-miRNA lentivirus?

While each anti-miRNA encodes a specific antisense RNA transcript, a negative control anti-miRNA (**a-miRNA-Neg-control**) lentivirus encodes a designed control sequence with minimal non-specific effects.

4. Why use GenTarget's anti-miRNA lentivirus:

GenTarget Inc

- The strongest and most stable anti-miRNA expression. Lentivirus can be effectively transduced into most dividing and non-dividing cells, and anti-miRNA can integrate into the host cell's genome for stable, long-term expression.
- **Compatibility.** The human H1 and human U6 promoters are active in almost all types of mammalian cells.
- **Constitutive or inducible expression.** GenTarget's lentivirus can be used for constitutive high shRNA expression, or optionally, for tetracycline inducible shRNA expression (with the H1 promoter only).
- **Full coverage.** You can order any anti-miRNAs listed in **miRBase** from any species; simply provide us the miRNA ID or the mature sequence, or your own validated anti-miRNA sequences.
- **Options for antibiotic markers.** Select from puromycin, blasticidin, or dual markers such as GFP-puromycin, GFP-blasticidin, RFP-puromycin, RFP-blasticidin.



Transduction Protocols:

1. Transduction Protocol for Adhesive cells:

Note: Pre-made lentivirus is provided ready to use, so it can be simply added into your cell culture; the amount of virus to add depends on cell type. For quick transduction, add 50 μ l of virus into each well of 24-well-plate where cell density is 50% to 75%. After 72 hours (no need to change medium), visualize positive transduction rate by fluorescence microscopy. For stable cell line generation, pass cells into medium containing antibiotic or perform fluorescence cell sorting followed by antibiotic selection.

Day 0:

Seed cells in complete medium at the appropriate density and incubate overnight.

Note: at the time of transduction, cells should be 50%-75% confluent. For example, seed HeLa cells at 0.5×10^{5} /ml x 0.5ml in a well of a 24-well plate.

Day 1:

- Thaw the pre-made lentiviral stock at room temperature and add the appropriate amount of virus stock to obtain the desired MOI.
- Return cells to 37°C, CO₂ incubator. Do nothing.
 Note: Try to avoid freezing and thawing. If you do not use all of the virus at one time, you may re-freeze the virus at -80 °C for future use; virus titer will decrease by ~10% for each freeze/thaw cycle.

Day 3:

At 48hr~72hr (Depend upon cell type) after transduction, check the transduction rate by fluorescence microscopy or calculate the exact transduction rate by flow cytometry (FACS or Guava).

Day 3 + (optional):

Sort transduced cells by FACS, or select by antibiotic killing. A pilot experiment should be done to determine the antibiotic's kill curve for your specific cell line (refer to the pertinent literature on generation of stable cell lines).

2. Transduction Protocol for Suspension Cells:



Grow cells in complete suspension culture medium; use a shaking flask in a CO^2 incubator if required.

Measure cell density (not grow over 3 million/ml), measured viability should be > 90%. Dilute cells into 1×10^6 cell/ml in complete medium.

Day 1:

• Thaw lentiviral particles at room temperature.

GenTarget Inc

- Add premade lentiviral particles into the diluted cells at a ratio of: 50 to 100 μ l virus per 0.5 ml of cells (Note: depending on cell type, you may need to use more or less virus).
- Grow cells in a shaking flask in a CO2 incubator.

Day 2:

At 24 hours after transduction, add an equal amount of fresh medium containing. Continue growing cells in CO2 incubator.

Day 3+:

At 48 hour to 72 hours (Depend upon cell type) after transduction, check fluorescence with a fluorescence microscope or calculate the transduction efficiency using a cell sorter such as FACS or Guava. Pass cells into 0.5 million/ml density in completed medium containing the corresponding antibiotic (**Note:** amount of antibiotic depends on cell type. A killing curve must pre-established). Sort for fluorescence positive cells and maintain antibiotic selection to generate a stable cell line.

Note: Filter wavelength settings: GFP filter: ~Ex450-490 ~Em525; RFP filter: ~Ex558 ~Em583;

Warranty:

This product is for research use only. It is warranted to meet its quality as described when used in accordance with its instructions. GenTarget disclaims any implied warranty of this product for particular application. In no event shall GenTarget be liable for any incidental or consequential damages in connection with the products. GenTarget's sole remedy for breach of this warranty should be, at GenTarget's option, to replace the products.



Safety Precaution:

Gentarget lentiviral particles adapts must advanced lentiviral safety features (using the third generation vectors with self-inactivation SIN-3UTR), and the premade lentivirus is replication incompetent. However, please use extra caution when using lentiviral particles. Use the lentiviral particles in Biosafety II cabinet. Wear glove all the time when handling Lentiviral particles! Please refer CDC and NIH's guidelines for more details regarding to safety issues.

References:

- 1. Lee RC, Feinbaum RL, Ambros V (December 1993). "The C. elegans heterochronic gene lin-4 encodes small RNAs with antisense complementarity to lin-14". Cell 75 (5): 843–54;
- 2. Lewis BP, Burge CB, Bartel DP (2005). "Conserved seed pairing, often flanked by adenosines, indicates that thousands of human genes are microRNA targets". Cell 120 (1): 15–20;
- 3. Hammond, S. M. Dicing and slicing: The core machinery of the RNA interference pathway. FEBS Lett. 2005, 579, 5822–5829.;
- 4. NIH Guidelines for **Biosafety Considerations for Research with Lentiviral Vectors**. (Link).

Related Products:	GenTarget's	pre-made	lentivirus	product	categories.
--------------------------	-------------	----------	------------	---------	-------------

Product Category	Product Description (please click into each category's page)
<u>Pathway</u> <u>Reporter</u>	Repoter Lentivirus for all kinds of pathway screening assays
<u>Cell</u> Immortalization	Lentivirus for cell immortalization: Large T-antigen, hTERT, EBNA1/EBNA2, HpV16-E6/E7, Adenovial E1A, Kras_G12V, HOXA9, et al.
<u>ImmunoOncology</u> <u>Research</u>	Lentivirus products for immuno therapy research: CAR and TCR; Assay Cell Lines for T-cell targeted killing assay and other cell-based assays; over-expression lentivirus products for the immune response targets; Cell surface antigens (CDs); immune checkpoint / Receptors; CRISPR gene Repair and knock-IN lentivirus; CRISPR knockout lentivirus;
<u>CAR-T, TCR</u> <u>Lentivirus</u>	CARs Lentivirus: Anti-CD19 /CD20 /CD22 /BCMA /hHER2 /HLA-A2 /TGFβ; TCRs : MART-1/ NY-ESO1/ CD1d-α-GalCer/ TRαV3-F2A-TRβV5-6;
<u>CRISPR Gene</u> <u>Editing</u>	Preamde lentivirus express humanzied wild-type Cas9 endonuclease, the dCas9, gRNAs, CRISPR gene





Product	Product Description	
Category	(please click into each category's page)	
	editing research	
Epigenomic:	"dCas9-Protein" fusion Lentivirus for epigenomic	
CRISPRI and	modification, resulted in CRISPR interference (CRISPRI)	
CRISPRA	or activation (CRISPRA).	
Cell-Specific	or fluorescent reporter (firefly Luciferase Renilla	
Reporter	luciferase. RFP or GFP fluorescent marker) under a	
	tissue specific promoter	
Infectious	Llentivirus that express all kinds of infectious antigens	
<u>Antigens</u>	with C-term 6His-tag.	
Virus Like	Lentiviral Like Particles, pseudo-typed with a different	
Particles (VLP)	envelope proteins.	
Non-integrating	Integration Defective Lentivirus, express different	
LV	targets for transient expression without the unwanted	
	insertional mutagenesis.	
<u>shRNA</u>	Knockdown verifeid and customized shRNA lentivirus for	
Knockdown	target knockdown,	
microRNA	Premade lentivirus expression human or mouse	
lentivirus	precursor miRNA. And anti-miRNA lentivector and	
Anti miNA	VIrus for numan and mouse mikina.	
Anu-mina Jentivirus		
Human and	Premade lentivirus expressin a human, mouse or rat	
<u>mouse ORFs</u>	gene with RFP-Blasticicin fusion dual markers.	
Luciferase	Premade lentivirus for all kinds of luciferase protein	
<u>expression</u>	expression: firefly and Renilla, Red-Luc and more,	
Elugroscopt	With different antibiotic selection markers.	
Markers	proteins: GED_RED_CED_RED_VED_piRED_upstable_GED	
<u>markers</u>	and others.	
Luminescent	Lentivirus express Nano-Latern as Bio-probes for in vivo	
Imaging	imaging of sub-cellular structural organization and	
	dynamic processes in living cells and organisms	
Sub-cellular	Lentivirus contain a well-defined organelle targeting	
<u>Imaging</u>	signal fusioned to a fluorescent protein, great tools for	



GenTarget Inc

Product	Product Description	
Category	(please click into each category's page)	
	live-cell imaging and for dynamic investigation of sub-	
	cellular signal pathways.	
<u>Cytoskeleton</u>	A fluorescent marker (GFP, RFP or CFP) fusion with a	
<u>Imaging</u>	cellular structure protein, provides a convenient tool	
	for visualization of cytoskeletal structure	
Unstable GFP	Lentivirus express the the destabilized GFP (uGFP)	
	which provides fast turnover responses in signal	
	pathway assay and in knockdown / knockout detection	
near-infrared RFP	The near-infrared Red fluorescent (niRFP) expression	
	Lentiviurs provides the whole-body images with better	
Eluprocept ODE	Contrast and Drighter integes	
fusion	fusion target	
	Promode lentivirus for expressing nuclear normeant	
CPE recombinase	CPE recombinase with different flurescent and antibiotic	
	markers	
CRF Fln	Lentivirus expressing "LoxP-GEP-Stop-LoxP-REP" or	
ColorSwtich	"FRT-GFP-Stop-FRT-RFP" cassette, used to monitor the	
	CRE or Flp recombination event in vivo.	
	lentivirus expressing SEAP under different promoters	
SEAP Reporter	(TetCMV, EF1a, CAG, Ubc, mPGK, Actin-beta or a	
	signal pathway responsive promoter),	
	Premade lentivirus expressin TetR (tetracycline	
TetR Repressor	regulator) protein, the repressor protein for the	
	inducible expression system.	
	rt I A binds to the tetracycline operator element (TetO)	
<u>rtra expression</u>	OFF inducible system	
	Premde lentivirus for human and mouse iPS (Myc	
iPS factors	NANOG OCTA SOX2 FLF4) factors with different	
	fluorescent and antibitoic markers	
LacZ expression	Express different full length B- galactosidase	
	(lacZ) with different selection markers	
	Premade negative control lentivirus with different	
Negative control	markers: serves as the negative control of lentivurs	
lentiviruses	treatment, for validation of the specificity of any	
	lentivirus target expression effects.	
Other Enzyme	Ready-to-use lentivirus, expressing a specific enzymes	





Product Category	Product Description (please click into each category's page)
expression	with different selection markers.
<u>Ultra titer</u> lentivirus	Ultra-titer lentivirus used for the hard-to-transduced cells and for in vivo manipulation of sperm cells, or stem cells.